AvaLight-DHc Full-range Compact Light Source

AvaLight-DHc



Get the best out of two worlds with the AvaLight-DHc. It has both a deuterium light source and a halogen light source, providing you with adequate light between 200 and 2500 nm for nearly all absorbance chemistry applications. Deuterium emits light between 200 and 550 nm, where the halogen takes over up to 2500 nm. Coupling this lightsource to the rest of your spectroscopy system is easy with the SMA connector.

This light source is recommended in settings with large fiber cables or directattachment to a cuvette holder such as the CUV-DA, due to its relatively low output energy. The integrated TTL-shutter makes saving a dark measurement very simple in combination with AvaSoft (extra IC-DB26-2 needed). Optionally the AvaLight-DHc is available in a rack-mountable version, to be used in the 19" rack or the 9.5" desktop system.

- Combined Deuterium-Halogen
- Integrated TTL-shutter

A direct-attach cuvette holder CUV-DA (see section accessories) is available for fluorescence or absorbance measurements.

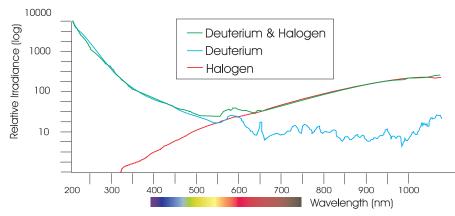


Figure 8 Spectral output of AvaLight-DHc

Technical Data

	Deuterium Light Source		Halogen Light Source
Wavelength Range	200 - 400 nm		400 - 2500 nm
Stability	< 1 mAU		< 1 mAU
Warm-up time	8 min		1 min
Drift	< 0.25% / h		< 0.25% / h
Optical Power in 600 µm fiber	0.2 μWatt		7 μWatt
Lamp Lifetime	1000 hours		2000 hours
Temp. Range	5°C - 35°C		
Power Supply	12VDC / 450 mA		
Dimensions, weight	175 x 110 x 44 mm, 570 g		
Lifetime shutter	1.000.000 cycles (typical)		




nordics@amstechnologies.com azpect.amstechnologies.com





Ordering Information

AvaLight-DHc	Compact Deuterium Halogen Light Source with TTL Shutter	
IC-DB26-2	• Interface cable AvaSpec-USB2/EVO platform to AvaLight-DHc-TTL-shutter	
AvaLight-DHc-RM	• Rackmount Compact Deuterium Halogen Light Source with TTL Shutter	
AvaLight-DHc-B	Compact Deuterium Halogen Replacement Bulb	
CUV-DA	• Direct-attach cuvette holder for AvaLight-DHc/XE/LED	
PS-12V/1.0A	• Power supply 100-240VAC/12VDC, 1.0A for AvaLight-DHc	

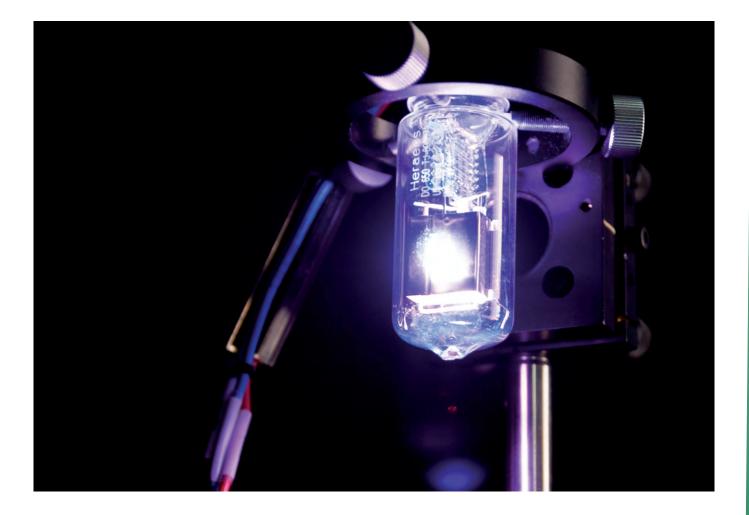

Three years limited warranty on all Avantes spectrometers, light sources and accessories





nordics@amstechnologies.com azpect.amstechnologies.com

Contact us

R